



Photo-reactive Crosslinkers

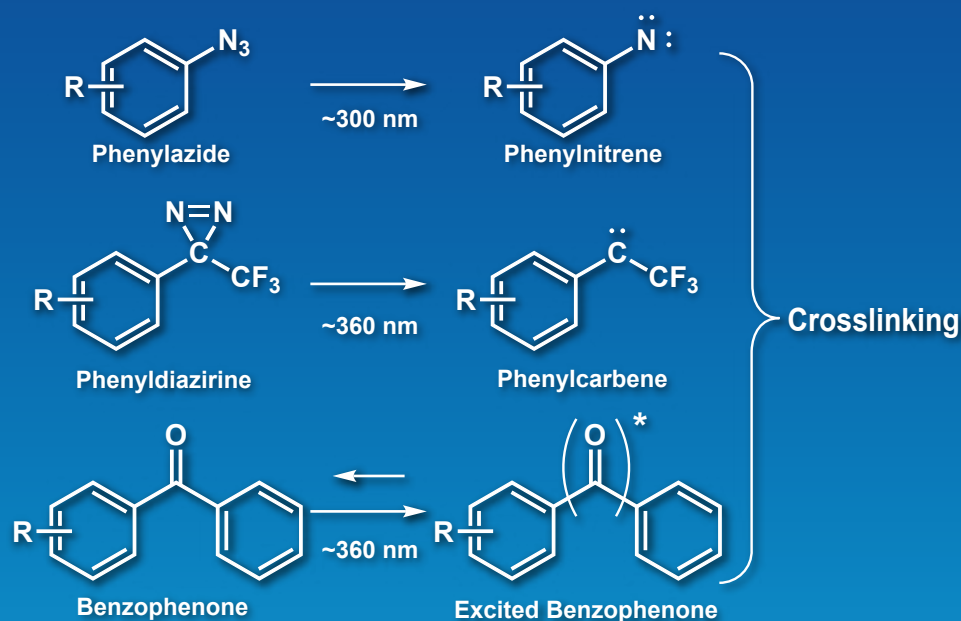

Some functional groups can be induced to react with target molecules by exposure to UV light. Photoreactive groups such as azides and diazirines are readily decomposed by UV light irradiation, resulting in highly reactive intermediates and the formation of covalent bonds with neighboring molecules. Within the context of molecular biology, photoaffinity labeling has prolific use in preparing ligands with photo-reactive units enabling the direct probing of target proteins.

Phenyldiazirines

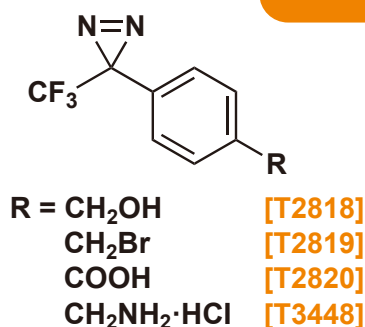

Phenyldiazirine generates a carbene unit by UV irradiation (<360 nm). Phenylcarbene can crosslink by short-time irradiation due to higher reactivity than nitrenes. Phenylcarbene is inactivated by water when neighboring target molecules are absent, and thus does not lead to non-specific crosslinking.

Applications

Reviews

- T. Tomohiro, M. Hashimoto, Y. Hatanaka, *Chem. Record* **2005**, 5, 385.
 M. Hashimoto, Y. Hatanaka, *Eur. J. Org. Chem.* **2008**, 2513.

Photoaffinity labeling

- Y. Kashiwayama, T. Tomohiro, K. Narita, M. Suzumura, T. Glumoff, J. K. Hiltunen, P. P. V. Veldhoven, Y. Hatanaka, T. Imanaka, *J. Biol. Chem.* **2010**, 285, 26315.
 E. W. S. Chan, S. Chattopadhyaya, R. C. Panicker, X. Huang, S. Q. Yao, *J. Am. Chem. Soc.* **2004**, 126, 14435.
 K. Matsuda, M. Ihara, K. Nishimura, D. B. Sattelle, K. Komai, *Biosci. Biotechnol. Biochem.* **2001**, 65, 1534.
 M. Wiegand, T. K. Lindhorst, *Eur. J. Org. Chem.* **2006**, 4841.

Photoaffinity microarray


- D. M. Dankbar, G. Gauglitz, *Anal. Bioanal. Chem.* **2006**, 386, 1967.
 S. Wei, J. Wang, D.-J. Guo, Y.-Q. Chen, S.-J. Xiao, *Chem. Lett.* **2006**, 35, 1172.
 N. Kanoh, S. Kumashiro, S. Simizu, Y. Kondoh, S. Hatakeyama, H. Tashiro, H. Osada, *Angew. Chem. Int. Ed.* **2003**, 42, 5584.

4-[3-(Trifluoromethyl)-3H-diazirin-3-yl]benzyl Alcohol
 4-[3-(Trifluoromethyl)-3H-diazirin-3-yl]benzyl Bromide
 4-[3-(Trifluoromethyl)-3H-diazirin-3-yl]benzoic Acid
 4-[3-(Trifluoromethyl)-3H-diazirin-3-yl]benzylamine Hydrochloride

200mg / 1g [T2818]
 200mg / 1g [T2819]
 200mg / 1g [T2820]
 200mg / 1g [T3448]

Phenylazides

Phenylazide generates a nitrene by UV irradiation (<300 nm). It is noted that azido groups tend to have less harmful effect on target analyte. Activation of the nitrene requires a shorter wavelength of UV light, and potential protein denaturation during long-period irradiation should be taken into consideration.

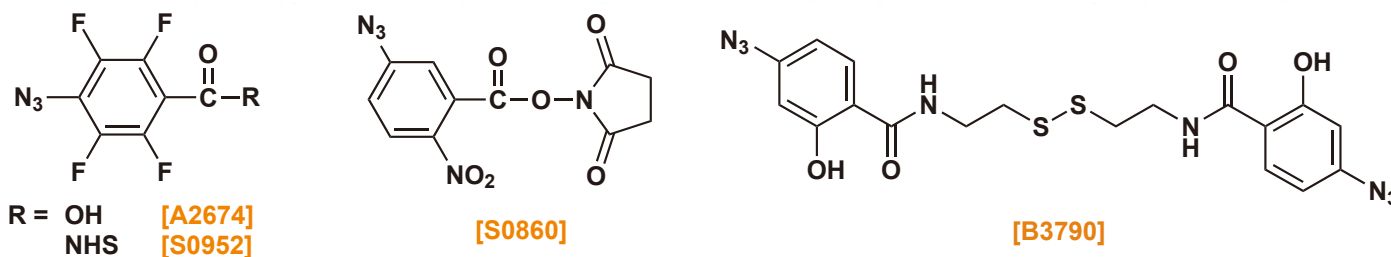

Applications

Photoaffinity labeling

J. F. W. Keana, S. Xiong Cai, *J. Org. Chem.* **1990**, 55, 3640.

Photoaffinity microarray

M. Thust, M. J. Schöning, J. Vetter, P. Kordos, H. Lüth, *Anal. Chim. Acta* **1996**, 323, 115.

Photo-crosslinking of protein complex

J. Rappsilber, S. Siniosoglou, E. C. Hurt, M. Mann, *Anal. Chem.* **2000**, 72, 267.

Photo-modification of carbon nanotube surface

S. J. Pastine, D. Okawa, B. Kessler, M. Rolandi, M. Llorente, A. Zettl, J. M. J. Fréchet, *J. Am. Chem. Soc.* **2008**, 130, 4238.

4-Azido-2,3,5,6-tetrafluorobenzoic Acid

1g [A2674]

4-Azido-2,3,5,6-tetrafluorobenzoic Acid N-Succinimidyl Ester

200mg / 1g [S0952]


5-Azido-2-nitrobenzoic Acid N-Succinimidyl Ester

10mg [S0860]

Bis[2-(4-azidosalicylamido)ethyl] Disulfide

10mg [B3790]

Benzophenone

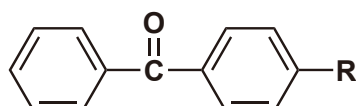

R = CO-NHS [S0863]

R = maleimide [M3259]

Benzophenone excited by UV irradiation (near 360 nm) to induce hydrogen abstraction from target molecules. The reaction efficiency remains high despite this due to the reverseability of the excited state. Additionally, photoexcited benzophenone is not water-reactive.

Applications

Photoaffinity labeling

G. F. Ross, P. M. Smith, A. McGregor, D. M. Turnbull, R. N. Lightowlers, *Bioconjugate Chem.* **2003**, 14, 962.

Y. Jung, J. M. Lee, J.-w. Kim, J. Yoon, H. Cho, B. H. Chung, *Anal. Chem.* **2009**, 81, 936.

Photoaffinity microarray

A. J. Hughes, A. E. Herr, *Proc. Natl. Acad. Sci. USA*, **2012**, 109, 21450.

4-Benzoylbenzoic Acid N-Succinimidyl Ester

200mg / 1g [S0863]

4-(N-Maleimido)benzophenone

50mg / 250mg [M3259]

For further information please refer to our website at www.TCIchemicals.com.

crosslinker



Ordering and Customer Service

TCI AMERICA

Tel : 800-423-8616 / 503-283-1681
Fax : 888-520-1075 / 503-283-1987
E-mail : Sales-US@TCIchemicals.com

TCI EUROPE N.V.

Tel : +32 (0)3 735 07 00
Fax : +32 (0)3 735 07 01
E-mail : Sales-EU@TCIchemicals.com

TCI Deutschland GmbH

Tel : +49 (0)6196 64053-00
Fax : +49 (0)6196 64053-01
E-mail : Sales-DE@TCIchemicals.com

Tokyo Chemical Industry UK Ltd.

Tel : +44 (0)1865 784560
Fax : +44 (0)1865 784561
E-mail : Sales-UK@TCIchemicals.com

TCI Chemicals (India) Pvt. Ltd.

Tel : 1800 425 7889 / 044-2262 0909
Fax : 044-2262 8902
E-mail : Sales-IN@TCIchemicals.com

梯希爱(上海)化成工业发展有限公司

Tel : 800-988-0390 / 021-67121386
Fax : 021-6712-1385
E-mail : Sales-CN@TCIchemicals.com

TOKYO CHEMICAL INDUSTRY CO., LTD.

Tel : +81 (0)3-5640-8878
Fax : +81 (0)3-5640-8902
E-mail : globalbusiness@TCIchemicals.com

Availability, price or specification of the listed products are subject to change without prior notice. Reproduction forbidden without the prior written consent of Tokyo Chemical Industry Co., Ltd.