



Glycosynthase

(Endo-M-N175Q)

An enzyme that adds whole sugar chains without breaking down products.


Increases efficiency of the sugar metastasis reaction


Applicable to Biosimilars

Making of a new functional sugar complex




Glycosynthase (Endo-M-N175Q) is an enzyme developed by Yamamoto, Umekawa, *et al.* through site-directed mutation of areas around the active center of Endo-M¹⁾ which is already marketed. Since the feature of Glycosynthase is efficient transglycosylation activity by using oxazoline derivatives as glycosyl donors while suppressing sugar hydrolysis activity, the resulting glycosylated products are obtained in high yield with less digestion of the products by the enzyme. Due to this feature Glycosynthase is expected to be applied as a useful tool in glycotechnology.

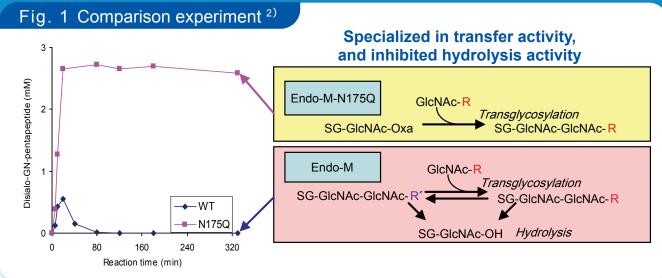

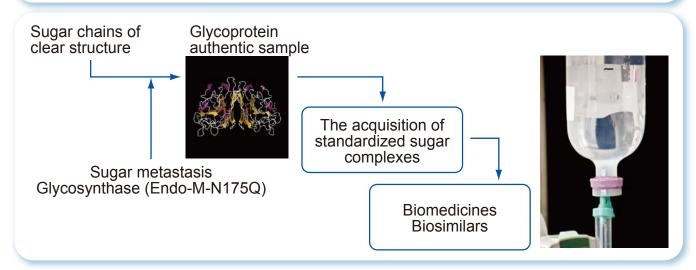


Umekawa and her colleagues caused transglycosylation reactions at the GlcNAc site of sperm antigen CD52 using oxazoline derivatives of the high-mannose type sugar chains or the complex type sugar chains as glycosyl donors²). They succeeded in obtaining glycosylated products in high yield of 84% and 76%, respectively. Moreover, they also achieved transglycosylation reactions using two biologically active blood-pressure-lowing peptides, PAMP12 and Substance P, as glycosyl acceptors and an oxazoline derivative of a complex type sugar chain containing sialic acids as a glycosyl donor in 95% and 98% yield, respectively³).

Fig. 2 Experiment example of the transglycosylation

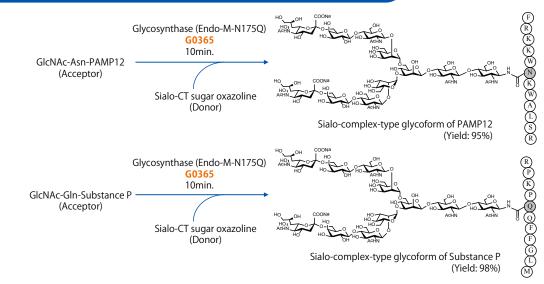

HPLC monitoring of the sugar metastasis reaction

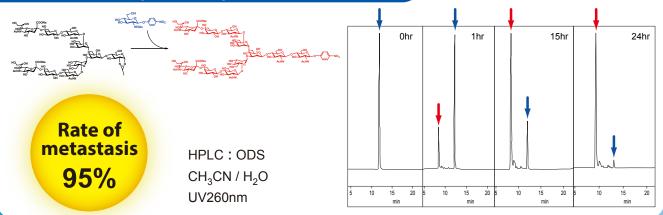

Realizing efficient transglycosylation reactions would also be useful for expansion into glycoprotein synthesis, such as the area of biosimilars.

Creation of new functional sugar complexes can be expected also.

Glycosynthase (Endo-M-N175Q)

from Mucor hiemalis expressed in Escherichia coli

Endo-B-N-Acotylalusosaminidaso (Endo-M

Endo-β-N-Acetylglucosaminidase (Endo-M)

from Mucor hiemalis expressed in Candida boidinii

Endo-M-W251N Recombinant

from Mucor hiemalis expressed in Escherichia coli

Sialylglycopeptide (SGP) Disialyloctasaccharide

100mU/vial [G0365]

100mU/vial [A1651]

500mU/vial [E1339]

10mg [S0523]

10mg [D4065]

- 1) K. Yamamoto, S. Kadowaki, J. Watanabe, H. Kumagai, Biochem. Biophys. Res. Commun. 1994, 203, 244.
- 2) M. Umekawa, T. Higashiyama, T. Tanaka, M. Noguchi, A. Kobayashi, S. Shoda, W. Huang, L-X. Wang, H. Ashida, K. Yamamoto, Biochim. Biophys. Acta, Gen. Subj. 2010, 1800, 1203.
- 3) M. Umekawa, C. Li, T. Higashiyama, W. Huang, H. Ashida, K. Yamamoto, L-X. Wang, J. Biol. Chem. 2010, 285, 511.

Ordering and Customer Service

TCI AMERICA

Tel : 800-423-8616 / 503-283-1681 Fax : 888-520-1075 / 503-283-1987 E-mail : Sales-US@TCIchemicals.com

TCI EUROPE N.V.

Tel : +32 (0)3 735 07 00 Fax : +32 (0)3 735 07 01 E-mail : Sales-EU@TCIchemicals.com

TCI Deutschland GmbH

Tel : +49 (0)6196 64053-00 Fax : +49 (0)6196 64053-01 E-mail : Sales-DE@TCIchemicals.com

Tokyo Chemical Industry UK Ltd.

Tel :+44 (0)1865 784560 Fax :+44 (0)1865 784561 E-mail : Sales-UK@TCIchemicals.com

E-mail: Sales-IN@TCIchemicals.com

TCI Chemicals (India) Pvt. Ltd.

Tel : 1800 425 7889 / 044-2262 0909 Fax : 044-2262 8902

E-mail: Sales-CN@TCIchemicals.com

:800-988-0390 / 021-67121386

梯希爱(上海)化成工业发展有限公司

Tel

TOKYO CHEMICAL INDUSTRY CO., LTD.
Tel : +81 (0)3-5640-8878

Fax : +81 (0)3-5640-8902 E-mail : globalbusiness@TCIchemicals.com

: 021-6712-1385

 $Availability, price \ or \ specification \ of \ the \ listed \ products \ are \ subject \ to \ change \ without \ prior \ notice. \ Reproduction \ forbidden \ without \ the \ prior \ written \ consent \ of \ Tokyo \ Chemical \ Industry \ Co., \ Ltd.$