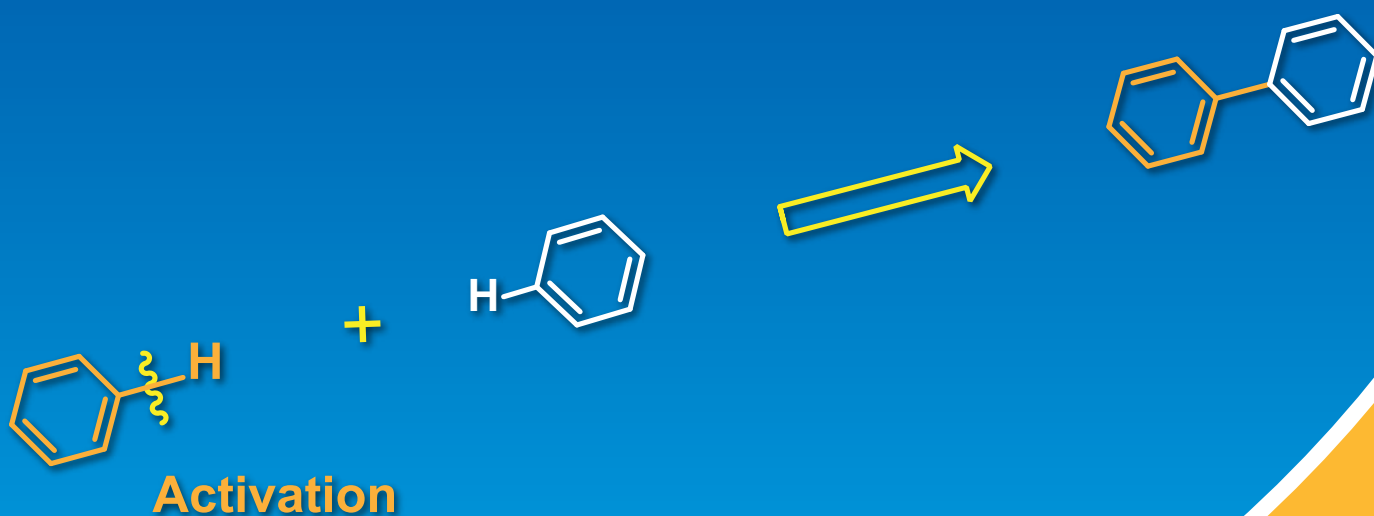


C-H Bond Activation Reaction



Metal Catalysts

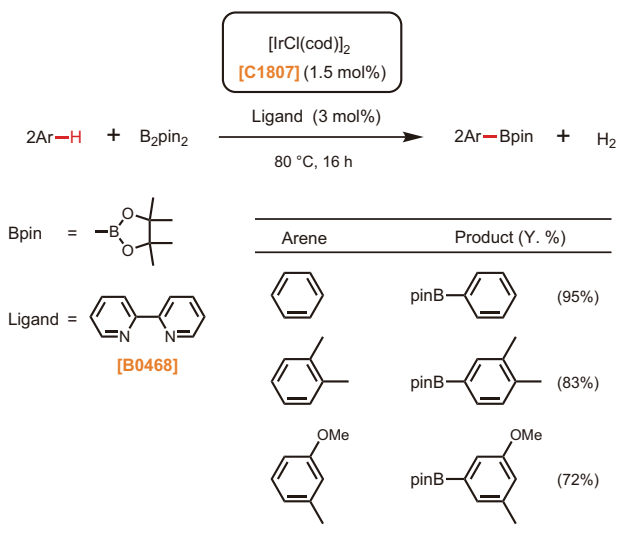
Ligands

Directing Group Introducing Agents

Additives

● Ir(I) Catalyst

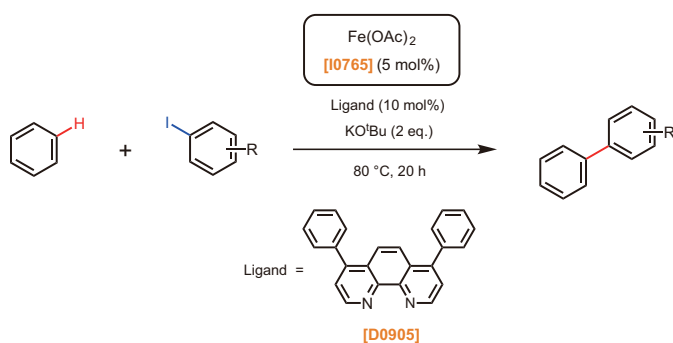
Miyaura, Ishiyama and Hargwig *et al.* have reported the direct C-H borylation in 2002.⁵⁾ This reaction is the most famous and practical example of C-H bond activation using iridium catalysts. Aryl borates had been synthesized by the reaction of aryl lithium or magnesium reagents with trialkyl borates so far, however, their method allowed a one-step preparation of alkyl borates in a simple manner.

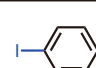
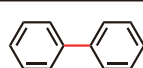
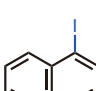
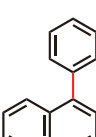
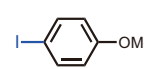
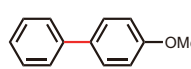
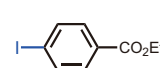
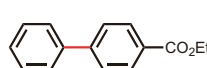
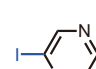
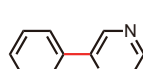


● Fe(II) Catalyst

Including palladium catalysts, which are frequently used for the Suzuki-Miyaura coupling reaction, transition metal catalysts, such as nickel or platinum, have been widely used for organic synthesis. However, the percentages of these metals in the earth's crust are extremely small, and their prices are rather expensive.⁶⁾ On the other hand, iron is abundant and less expensive, and therefore, more and more chemists have focused their attention to organic synthesis using iron compounds as a catalyst. Cross coupling reactions using iron catalysts have been reported.⁷⁾

For an example of C-H activation using iron catalysts, Charette *et al.* have reported the direct coupling reaction of benzene with aryl iodides using iron(II) acetate [I0765].⁸⁾ This reaction is highly cost-effective and environmentally friendly in the sense of using an iron catalyst, which is less expensive, and therefore, further development and applications are expected from the point of green chemistry.



Aryl Iodide	Product (Y. %)
	 (89%)
	 (60%)
	 (93%)
	 (40%)
	 (79%)

Thus, C-H bond activation has been widely studied as a new methodology of carbon-carbon and carbon-heteroatom bond formations, following a cross coupling reaction and olefin metathesis.

TCI offers a variety of transition metal catalysts, ligands, and activating reagents readily available for C-H bond activation as below.

References

- 1) S. Murai, F. Kakiuchi, S. Sekine, Y. Tanaka, A. Kamatani, M. Sonoda, N. Chatani, *Nature* **1993**, 366, 529.
- 2) K. L. Hull, M. S. Sanford, *J. Am. Chem. Soc.* **2007**, 129, 11904.
- 3) M. S. Chen, N. Prabakaran, N. A. Labenz, M. C. White, *J. Am. Chem. Soc.* **2005**, 127, 6970.
- 4) K. J. Fraunhofer, N. Prabakaran, L. E. Sirois, M. C. White, *J. Am. Chem. Soc.* **2006**, 128, 9032.
- 5) T. Ishiyama, J. Takagi, K. Ishida, N. Miyaura, N. R. Anastasi, J. F. Hartwig, *J. Am. Chem. Soc.* **2002**, 124, 390.
- 6) J. Emsley, in *The Elements*, 3rd ed., Oxford Univ. Press, New York, **1998**.
- 7) T. Nagano, T. Hayashi, *Org. Lett.* **2004**, 6, 1297; K. G. Dongo, H. Koh, M. Sau, C. L. L. Chai, *Adv. Synth. Catal.* **2007**, 349, 1015; T. Hatakeyama, M. Nakamura, *J. Am. Chem. Soc.* **2007**, 129, 9844.
- 8) F. Vallée, J. J. Mousseau, A. B. Charette, *J. Am. Chem. Soc.* **2010**, 132, 1514,

Metal Catalysts

Palladium Catalysts

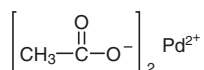
P1489 1g 5g

Palladium(II) Chloride
CAS RN: 7647-10-1

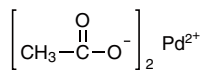
P1937 2g

Palladium(II) Bromide
CAS RN: 13444-94-5

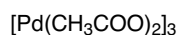
A1424 1g 5g

Palladium(II) Acetate
CAS RN: 3375-31-3

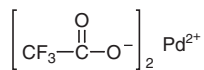
P2161 1g

Palladium(II) Acetate (Purified)
CAS RN: 3375-31-3

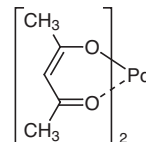
P2106 1g 5g

Palladium(II) Acetate Trimer
CAS RN: 53189-26-7

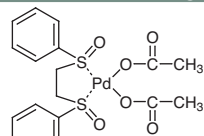
P1870 1g 5g

Palladium(II) Trifluoroacetate
CAS RN: 42196-31-6

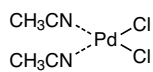
B2018 1g 5g

Palladium(II) Acetylacetonate
CAS RN: 14024-61-4

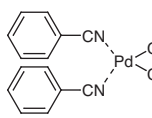
B3292 200mg 1g

White Catalyst
CAS RN: 858971-43-4

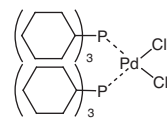
B1676 1g 5g

Bis(acetonitrile)palladium(II)
Dichloride
CAS RN: 14592-56-4

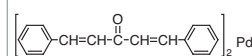
B1668 1g 5g

Bis(benzonitrile)-
palladium(II) Dichloride
CAS RN: 14220-64-5

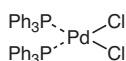
B2055 1g 5g

Bis(tricyclohexylphosphine)-
palladium(II) Dichloride
CAS RN: 29934-17-6

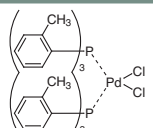
B1374 1g 5g

Bis(dibenzylideneacetone)-
palladium(0)
CAS RN: 32005-36-0

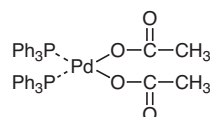
B1667 1g 5g 25g

Bis(triphenylphosphine)-
palladium(II) Dichloride
CAS RN: 13965-03-2

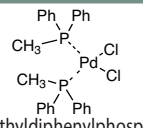
B2026 1g 5g

Bis(tri-*o*-tolylphosphine)-
palladium(II) Dichloride
CAS RN: 40691-33-6

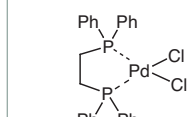
B2042 1g 5g

Bis(triphenylphosphine)-
palladium(II) Diacetate
CAS RN: 14588-08-0

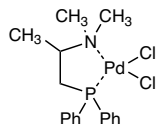
B2161 1g 5g

Bis(methyldiphenylphosphine)-
palladium(II) Dichloride
(*cis*- and *trans*-mixture)
CAS RN: 52611-08-2

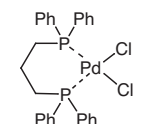
B2016 1g 5g

[1,2-Bis(diphenylphosphino)-
ethane]palladium(II) Dichloride
CAS RN: 19978-61-1

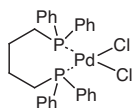
D5719 250mg

PdCl₂(alphos)
CAS RN: 85719-56-8

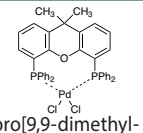
B2192 1g 5g

[1,3-Bis(diphenylphosphino)-
propane]palladium(II) Dichloride
CAS RN: 59831-02-6

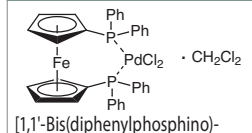
B2031 1g 5g

[1,4-Bis(diphenylphosphino)-
butane]palladium(II) Dichloride
CAS RN: 29964-62-3

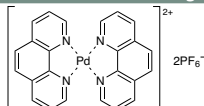
D4333 200mg 1g

Dichloro[9,9-dimethyl-
4,5-bis(diphenylphosphino)-
xanthene]palladium(II)
CAS RN: 205319-10-4

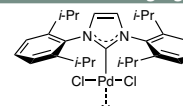
B2064 1g 5g 25g

[1,1'-Bis(diphenylphosphino)-
ferrocene]palladium(II) Dichloride
Dichloromethane Adduct
CAS RN: 95464-05-4

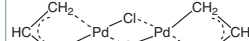
B5400 200mg 1g

Bis(1,10-phenanthroline)-
palladium(II)
Bis(hexafluorophosphate)
CAS RN: 113173-22-1

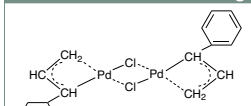
B6199 1g 5g 25g

Organ's Catalyst
CAS RN: 905459-27-0

A1479 500mg 1g

Allylpalladium(II) Chloride
Dimer
CAS RN: 12012-95-2

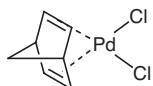
P2017 200mg

Palladium(II)(π -cinnamyl)-
Chloride Dimer
CAS RN: 12131-44-1

D2604 1g 5g

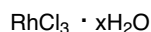
Dichloro(1,5-cyclooctadiene)-
palladium(II)
CAS RN: 12107-56-1

N0842 1g

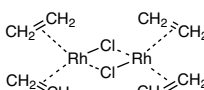
2,5-Norbornadiene
Palladium(II) Dichloride
CAS RN: 12317-46-3

Rhodium Catalysts

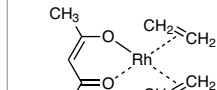
R0244 250mg

Rhodium(III) Chloride
Hydrate
CAS RN: 20765-98-4

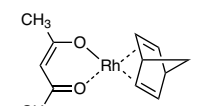
C2461 200mg

Chlorobis(ethylene)-
rhodium(I) Dimer
CAS RN: 12081-16-2

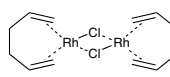
A2100 200mg

Acetylacetonatobis(ethylene)-
rhodium(I)
CAS RN: 12082-47-2

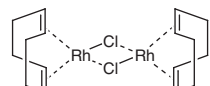
A2761 200mg 1g

(Acetylacetonato)-
(norbornadiene)rhodium(I)
CAS RN: 32354-50-0

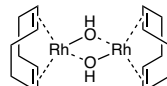
C3194 100mg

Chloro(1,5-hexadiene)-
rhodium(I) Dimer
CAS RN: 32965-49-4

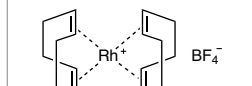
B1045 100mg 1g

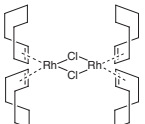
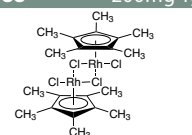
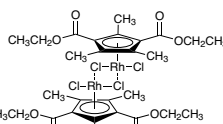
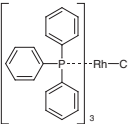
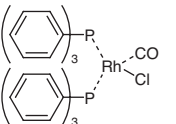
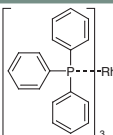
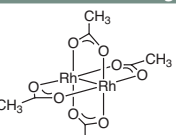
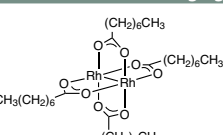
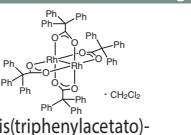
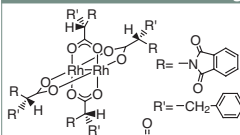
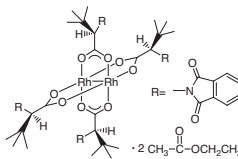
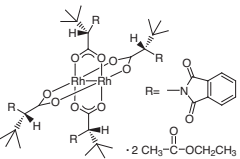
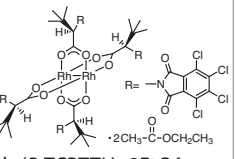
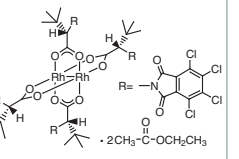
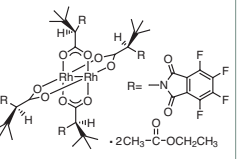
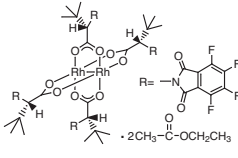
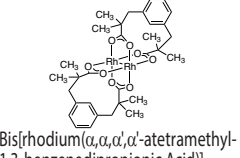
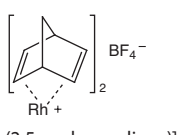
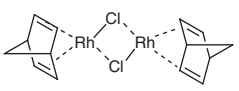
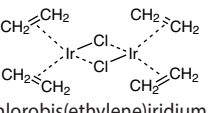
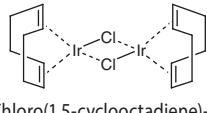
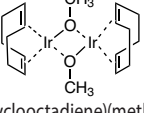
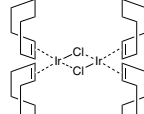
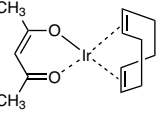
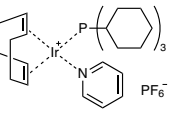
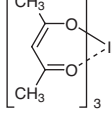
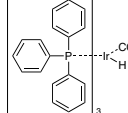
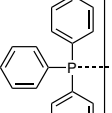
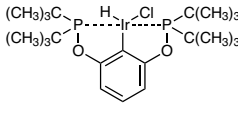
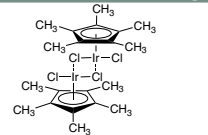
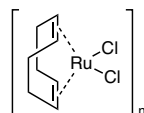
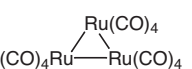
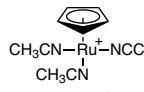
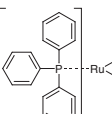
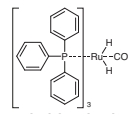
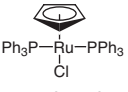
Chloro(1,5-cyclooctadiene)-
rhodium(I) Dimer
CAS RN: 12092-47-6

H1562 200mg 1g

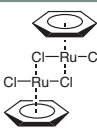
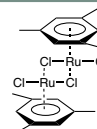
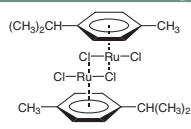
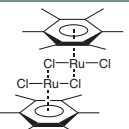
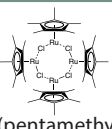
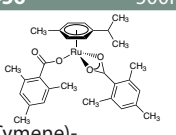
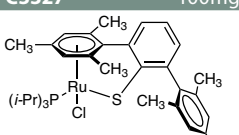
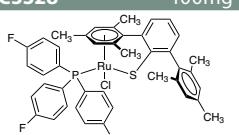
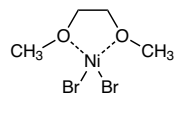
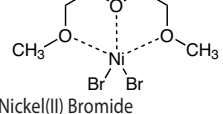
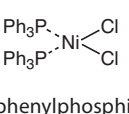
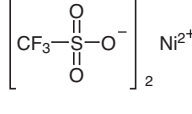
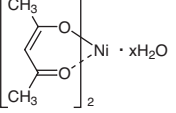
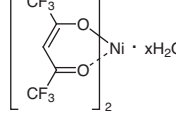
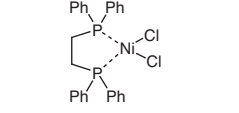
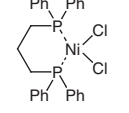
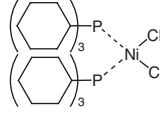
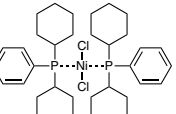
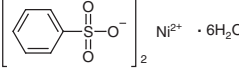
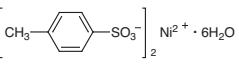
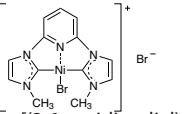
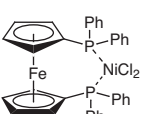
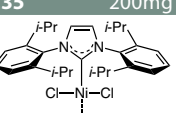
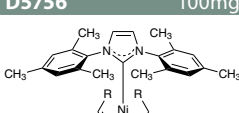
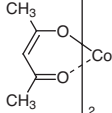
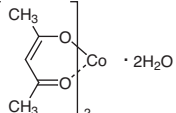
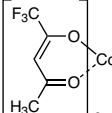
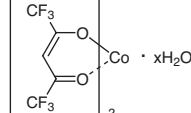
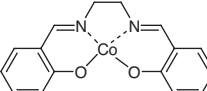
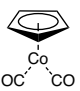
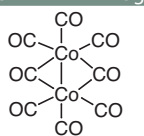
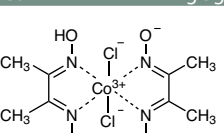
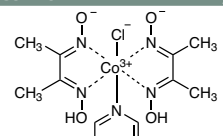
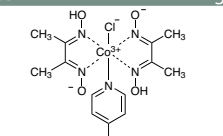
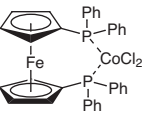
Hydroxy(1,5-cyclooctadiene)-
rhodium(I) Dimer
CAS RN: 73468-85-6

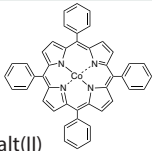
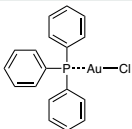
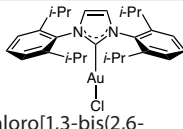
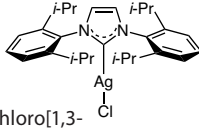
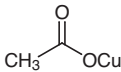
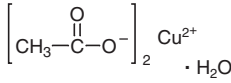
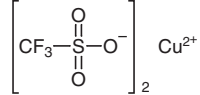
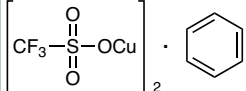
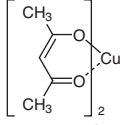
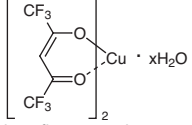
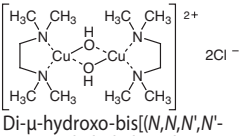
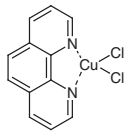
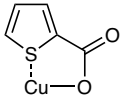
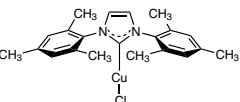
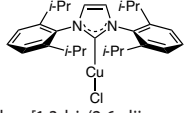
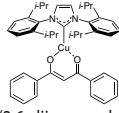
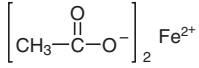
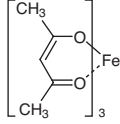
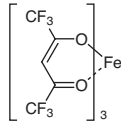
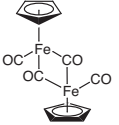
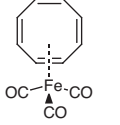
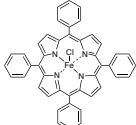
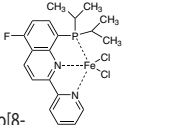
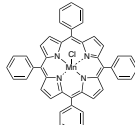
B3961 100mg 1g

Bis(1,5-cyclooctadiene)-
rhodium(I) Tetrafluoroborate
CAS RN: 35138-22-8

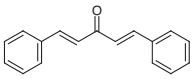
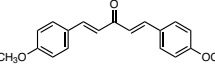
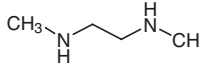


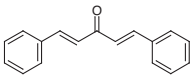
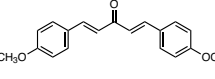
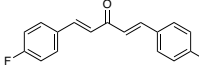
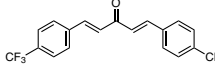
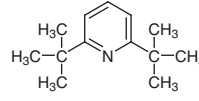
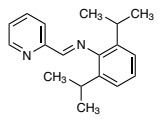
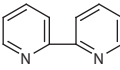
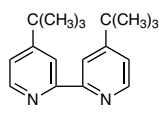
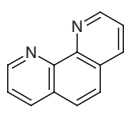
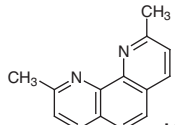
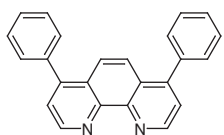
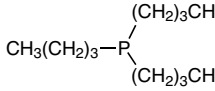
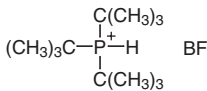
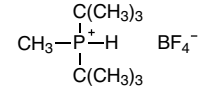
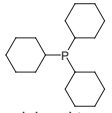
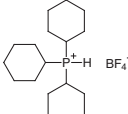
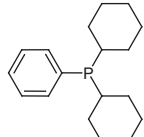
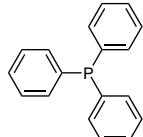
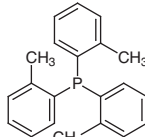
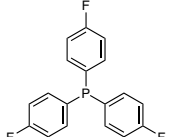
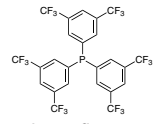
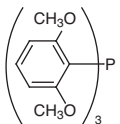
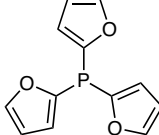
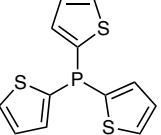
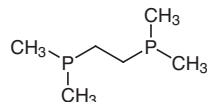
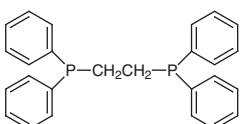
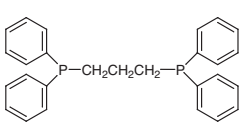
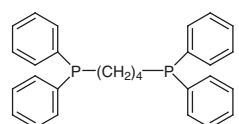
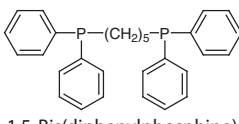
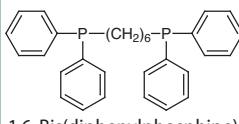
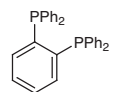
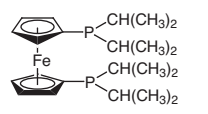
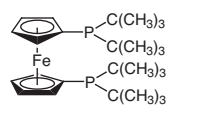
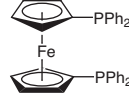
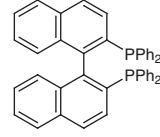
<p>C2253 100mg 500mg</p>  <p>Chlorobis(cyclooctene)rhodium(I) Dimer CAS RN: 12279-09-3</p>	<p>P1788 200mg 1g</p>  <p>(Pentamethylcyclopentadienyl)rhodium(III) Dichloride Dimer CAS RN: 12354-85-7</p>	<p>B6169 200mg</p>  <p>[Cp[*]RhCl₂]₂ CAS RN: 1352745-18-6</p>	<p>T0931 1g 5g</p>  <p>Tris(triphenylphosphine)rhodium(II) Chloride CAS RN: 14694-95-2</p>	<p>B1692 1g</p>  <p>Carbonylbis(triphenylphosphine)rhodium(II) Chloride CAS RN: 13938-94-8</p>
<p>C1383 1g 5g</p>  <p>Carbonylhydrido-tris(triphenylphosphine)rhodium(II) CAS RN: 17185-29-4</p>	<p>R0069 100mg 1g</p>  <p>Rhodium(II) Acetate Dimer CAS RN: 15956-28-2</p>	<p>R0161 200mg 1g</p>  <p>Rhodium(II) Octanoate Dimer CAS RN: 73482-96-9</p>	<p>T1544 100mg</p>  <p>Tetrakis(triphenylacetato)dirhodium(II) Dichloromethane Adduct CAS RN: 142214-04-8</p>	<p>T1551 100mg</p>  <p>Rh₂(S-PTPA)₄ EtOAc CAS RN: 131219-55-1</p>
<p>T2054 100mg</p>  <p>Rh₂(R-PTTL)₄ 2EtOAc</p>	<p>T2055 100mg</p>  <p>Rh₂(S-PTTL)₄ 2EtOAc</p>	<p>T2658 50mg</p>  <p>Rh₂(R-TCPTTL)₄ 2EtOAc CAS RN: 2001054-66-4</p>	<p>T2659 100mg</p>  <p>Rh₂(S-TCPTTL)₄ 2EtOAc</p>	<p>T2660 50mg</p>  <p>Rh₂(R-TFPTTL)₄ 2EtOAc</p>
<p>T2661 100mg</p>  <p>Rh₂(S-TFPTTL)₄ 2EtOAc</p>	<p>B4549 100mg 500mg</p>  <p>Bis[rhodium(α,α,α',α'-tetramethyl-1,3-benzenedipropionic Acid)] CAS RN: 819050-89-0</p>	<p>B2091 100mg</p>  <p>Bis[η-(2,5-norbornadiene)]rhodium(I) Tetrafluoroborate CAS RN: 36620-11-8</p>	<p>N0453 100mg</p>  <p>Norbornadiene Rhodium(I) Chloride Dimer CAS RN: 12257-42-0</p>	<p style="text-align: center;">Iridium Catalysts</p>
<p>I0616 1g 5g</p> <p style="text-align: center;">IrCl₃ · xH₂O</p> <p>Iridium(III) Chloride Hydrate CAS RN: 14996-61-3</p>	<p>C3041 100mg</p>  <p>Chlorobis(ethylene)iridium(I) Dimer CAS RN: 39722-81-1</p>	<p>C1807 250mg 1g</p>  <p>Chloro(1,5-cyclooctadiene)iridium(I) Dimer CAS RN: 12112-67-3</p>	<p>C2662 200mg 1g</p>  <p>(1,5-Cyclooctadiene)(methoxy)iridium(I) Dimer CAS RN: 12148-71-9</p>	<p>C2985 200mg</p>  <p>Chlorobis(cyclooctene)iridium(I) Dimer CAS RN: 12246-51-4</p>
<p>A2981 200mg 1g</p>  <p>(Acetylacetonato)(1,5-cyclooctadiene)iridium(I) CAS RN: 12154-84-6</p>	<p>C2824 100mg</p>  <p>Crabtree's Catalyst CAS RN: 64536-78-3</p>	<p>T2557 1g</p>  <p>Tris(2,4-pentanedionato)iridium(III) CAS RN: 15635-87-7</p>	<p>C3040 200mg 1g</p>  <p>Carbonylhydrido-tris(triphenylphosphine)iridium(II) CAS RN: 17250-25-8</p>	<p>C2252 200mg 1g</p>  <p>Vaska's Catalyst CAS RN: 14871-41-1</p>
<p>B5033 200mg</p>  <p>2,6-Bis(di-tert-butylphosphinoxy)phenylchlorohydroiridium(III) CAS RN: 671789-61-0</p>	<p>P1763 1g</p>  <p>(Pentamethylcyclopentadienyl)iridium(III) Dichloride Dimer CAS RN: 12354-84-6</p>	<p style="text-align: center;">Ruthenium Catalysts</p>	<p>R0074 1g 5g</p> <p style="text-align: center;">RuCl₃ · xH₂O</p> <p>Ruthenium(III) Chloride Hydrate CAS RN: 14898-67-0</p>	<p>D4792 1g 5g</p>  <p>[RuCl₂(COD)]_n CAS RN: 50982-12-2</p>
<p>T2181 100mg 1g</p>  <p>Triruthenium Dodecacarbonyl CAS RN: 15243-33-1</p>	<p>T3079 200mg</p>  <p>Tris(acetonitrile)cyclopentadienylruthenium(II) Hexafluorophosphate CAS RN: 80049-61-2</p>	<p>D1997 1g 5g</p>  <p>Tris(triphenylphosphine)ruthenium(II) Dichloride CAS RN: 15529-49-4</p>	<p>C2251 250mg 1g</p>  <p>Carbonyl(dihydrido)tris(triphenylphosphine)ruthenium(II) CAS RN: 25360-32-1</p>	<p>C2201 1g 5g</p>  <p>Cyclopentadienylbis(triphenylphosphine)ruthenium(II) Chloride CAS RN: 32993-05-8</p>

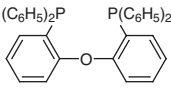
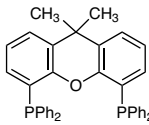
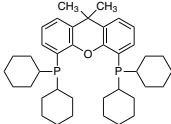
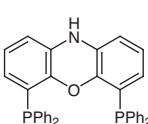
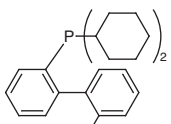
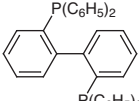
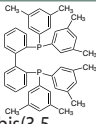
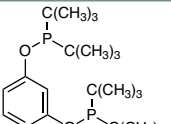
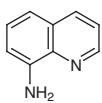
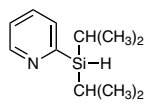
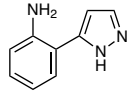
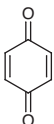
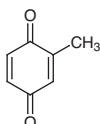
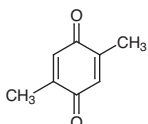
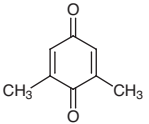
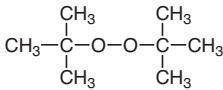
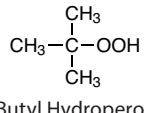
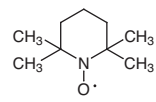
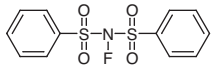
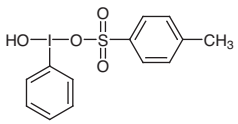
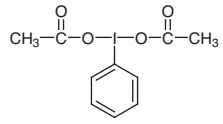
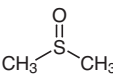
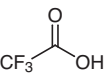
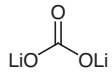
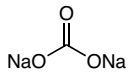
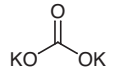
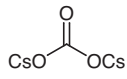
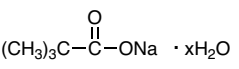
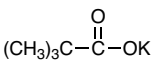
C-H Bond Activation Reaction

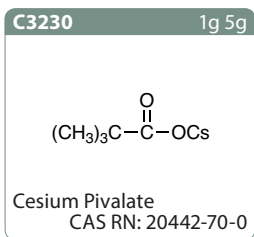
B1902 1g 5g  Benzeneruthenium(II) Chloride Dimer CAS RN: 37366-09-9	D5524 250mg 1g  Mesityleneruthenium(II) Chloride Dimer CAS RN: 52462-31-4	D2751 1g 5g  Dichloro(<i>p</i> -cymene)ruthenium(II) Dimer CAS RN: 52462-29-0	H1010 1g  (Hexamethylbenzene)-ruthenium(II) Dichloride Dimer CAS RN: 67421-02-7	C3042 200mg 1g  Chloro(pentamethylcyclopentadienyl)-ruthenium(II) Tetramer CAS RN: 113860-07-4	
C3456 500mg  (<i>p</i> -Cymene)-bis(mesitylcarboxylato)-ruthenium(II) CAS RN: 1251667-99-8	C3327 100mg  [(DmpSR)RuCl(P-(<i>i</i> -Pr) ₃)] CAS RN: 1621182-04-4	C3328 100mg  [(DmpSR)RuCl(P-(4-Fluorophenyl) ₃)] CAS RN: 1420299-84-8	<h2>Nickel Catalysts</h2>		N0850 25g 500g NiCl ₂ Nickel(II) Chloride Anhydrous CAS RN: 7718-54-9
N0851 25g 500g NiCl ₂ · 6H ₂ O Nickel(II) Chloride Hexahydrate CAS RN: 7791-20-0	N1050 1g 10g  NiBr (dme) CAS RN: 28923-39-9	N1049 1g 10g  Nickel(II) Bromide 2-Methoxyethyl Ether Complex CAS RN: 312696-09-6	B1571 10g 100g  Bis(triphenylphosphine)-nickel(II) Dichloride CAS RN: 14264-16-5	N0861 1g 5g  Nickel(II) Triflate CAS RN: 60871-84-3	
N0096 25g 100g 500g  Bis(2,4-pentanedionato)-nickel(II) Hydrate CAS RN: 120156-44-7	H0558 1g 5g  Bis(hexafluoroacetylacetonato)-nickel(II) Hydrate CAS RN: 14949-69-0	B2225 1g 5g 25g  [1,2-Bis(diphenylphosphino)ethane]nickel(II) Dichloride CAS RN: 14647-23-5	B1313 5g 25g  [1,3-Bis(diphenylphosphino)propane]dichloronickel(II) CAS RN: 15629-92-2	B3534 1g 5g  Bis(tricyclohexylphosphine)-nickel(II) Dichloride CAS RN: 19999-87-2	
D5369 250mg  Dichlorobis(dicyclohexylphenylphosphine)nickel(II) CAS RN: 19232-03-2	B0034 25g  Nickel(II) Benzenesulfonate Hexahydrate CAS RN: 39819-65-3	T0276 5g 25g  Nickel(II) <i>p</i> -Toluenesulfonate Hexahydrate CAS RN: 6944-05-4	B3354 1g  Bromo(2,6-pyridinediyl)-bis(3-methyl-1-imidazolyl-2-ylidene)nickel Bromide CAS RN: 894102-11-5	B2226 1g 5g  [1,1'-Bis(diphenylphosphino)ferrocene]nickel(II) Dichloride CAS RN: 67292-34-6	
B3235 200mg 1g  [1,3-Bis(2,6-diisopropylphenyl)-imidazol-2-ylidene]triphenylphosphine Nickel(II) Dichloride CAS RN: 903592-98-3	D5756 100mg  Ni(I)(Mes)(Di- <i>tert</i> -butyl Fumarate) ₂ CAS RN: 2230140-59-5	<h2>Cobalt Catalysts</h2>		C2388 250g CoCl ₂ · 6H ₂ O Cobalt(II) Chloride Hexahydrate CAS RN: 7791-13-1	B2681 25g  Acetylacetonate Cobalt(II) Salt CAS RN: 14024-48-7
C0373 25g 500g  Bis(2,4-pentanedionato)-cobalt(II) Dihydrate CAS RN: 123334-29-2	T0746 1g 5g  Bis(trifluoro-2,4-pentanedionato)cobalt(II) CAS RN: 16092-38-9	H0553 5g  Bis(hexafluoroacetylacetonato)-cobalt(II) Hydrate CAS RN: 19648-83-0	S0318 25g 100g 500g  Salcomine CAS RN: 14167-18-1	D4940 2g  Dicarbonyl-cyclopentadienylcobalt(I) CAS RN: 12078-25-0	
D3213 5g 25g  Dicobalt Octacarbonyl (stabilized with 1-5% Hexane) CAS RN: 10210-68-1	D5924 1g 5g  Co(dmgh) ₂ Cl ₂ CAS RN: 23638-66-6	C3718 1g  Co(dmgh) ₂ PyCl CAS RN: 23295-32-1	C3711 1g  Co(dmgh) ₂ (DMAP)Cl CAS RN: 483979-48-2	B3374 1g 5g  [1,1'-Bis(diphenylphosphino)ferrocene]cobalt(II) Dichloride CAS RN: 67292-36-8	

C3579 1g 5g  Cobalt(II) Tetraphenylporphyrin CAS RN: 14172-90-8	Gold Catalysts	T2994 200mg 1g  (Triphenylphosphine)- gold(I) Chloride CAS RN: 14243-64-2	C2405 200mg  Chloro[1,3-bis(2,6- diisopropylphenyl)- imidazol-2-ylidene]gold(I) CAS RN: 852445-83-1	Silver Catalysts
S0463 5g 25g AgSbF ₆ Silver Hexafluoroantimonate(V) CAS RN: 26042-64-8	S0981 1g 5g AgPF ₆ Silver Hexafluorophosphate CAS RN: 26042-63-7	S0898 1g 5g Ag ⁺ (CF ₃ SO ₂) ₂ N ⁻ Silver Triflimide CAS RN: 189114-61-2	C2373 200mg 1g  Chloro[1,3- bis(2,6-diisopropylphenyl)- imidazol-2-ylidene]silver CAS RN: 873297-19-9	Copper Catalysts
T2666 1g 5g 25g (CH ₃ CN) ₄ Cu ⁺ BF ₄ ⁻ Tetrakis(acetonitrile)copper(I) Tetrafluoroborate CAS RN: 15418-29-8	T2665 5g (CH ₃ CN) ₄ Cu ⁺ PF ₆ ⁻ Tetrakis(acetonitrile)copper(I) Hexafluorophosphate CAS RN: 64443-05-6	T3905 1g 5g (CH ₃ CN) ₄ Cu ⁺ CF ₃ SO ₃ ⁻ Tetrakis(acetonitrile)copper(I) Triflate CAS RN: 58452-28-1	C1952 25g 300g CuCN Copper(I) Cyanide CAS RN: 544-92-3	C2162 25g 100g 500g CuCl Copper(I) Chloride CAS RN: 7758-89-6
C2161 25g 100g 500g CuBr Copper(I) Bromide CAS RN: 7787-70-4	C2163 25g 100g 500g CuI Copper(I) Iodide CAS RN: 7681-65-4	A1540 5g 25g  Copper(I) Acetate CAS RN: 598-54-9	C2346 25g 500g  Copper(II) Acetate Monohydrate CAS RN: 6046-93-1	T1292 5g 25g  Copper(II) Trifluoromethanesulfonate CAS RN: 34946-82-2
T1442 1g 5g  Copper(I) Triflate Benzene Complex CAS RN: 42152-46-5	C0384 25g 250g  Bis(2,4-pentanedionato)- copper(II) CAS RN: 13395-16-9	H0554 1g 5g  Bis(hexafluoroacetylacetonato)- copper(II) Hydrate CAS RN: 14781-45-4	D2542 5g 25g  Di-μ-hydroxo-bis[(N,N,N',N'- tetramethylethylenediamine)- copper(II)] Chloride CAS RN: 30698-64-7	D3891 1g 5g  Dichloro(1,10- phenanthroline)copper(III) CAS RN: 14783-09-6
C2312 1g 5g  CuTC CAS RN: 68986-76-5	C2422 200mg 1g  Chloro(1,3-dimesitylimidazol- 2-ylidene)copper(I) CAS RN: 873779-78-3	C2304 200mg 1g  Chloro[1,3-bis(2,6-diisopropyl- phenyl)imidazol- 2-ylidene]copper(I) CAS RN: 578743-87-0	B3351 200mg 1g  [1,3-Bis(2,6-diisopropylphenyl)- imidazol-2-ylidene](1,3-diphenyl- 1,3-propanedionato)copper(I) CAS RN: 920739-11-3	Iron Catalysts
I0765 5g 25g  Iron(II) Acetate CAS RN: 3094-87-9	I0079 25g 100g 500g  Iron(III) Acetylacetonate CAS RN: 14024-18-1	H0555 1g  Tris(hexafluoroacetylacetonato)- iron(III) CAS RN: 17786-67-3	C1592 5g  Cyclopentadienyliron Dicarboxyl Dimer CAS RN: 12154-95-9	T1775 1g  Tricarbonyl(cyclooctatetraene)- iron CAS RN: 12093-05-9
I0937 200mg 1g  Iron(III) Tetraphenylporphyrin Chloride CAS RN: 16456-81-8	D5886 100mg  Dichloro[8- (diisopropylphosphino)-5-fluoro- 2-(2-pyridinyl)quinoline]iron(II) CAS RN: 2247605-87-2	Manganese Catalysts	M2621 1g 5g  Manganese(III) Tetraphenylporphyrin Chloride CAS RN: 32195-55-4	D5880 1g [Mn(CO) ₅] ₂ Manganese Carbonyl CAS RN: 10170-69-1

Ligands

<p>D0903 25g 250g</p>  <p><i>trans,trans</i>-Dibenzylideneacetone CAS RN: 35225-79-7</p>	<p>B4467 200mg 1g</p>  <p><i>trans,trans</i>-Bis(4-methoxybenzylidene)acetone CAS RN: 37951-12-5</p>	<p>D0720 5mL 25mL</p>  <p><i>N,N'</i>-Dimethylethylenediamine CAS RN: 110-70-3</p>	<p>N0166 25g 400g</p>  <p>2-Norbornene CAS RN: 498-66-8</p>	<p>N0346 25mL 100mL 500mL</p>  <p>2,5-Norbornadiene (stabilized with BHT) CAS RN: 121-46-0</p>
<p>D0903 25g 250g</p>  <p><i>trans,trans</i>-Dibenzylideneacetone CAS RN: 35225-79-7</p>	<p>B4467 200mg 1g</p>  <p><i>trans,trans</i>-Bis(4-methoxybenzylidene)acetone CAS RN: 37951-12-5</p>	<p>B2283 5g 25g</p>  <p><i>trans,trans</i>-Bis(4-fluorobenzylidene)acetone CAS RN: 53369-00-9</p>	<p>B4468 200mg 1g</p>  <p><i>trans,trans</i>-Bis[4-(trifluoromethyl)benzylidene]acetone CAS RN: 103836-71-1</p>	<p>D1804 5g 25g</p>  <p>2,6-Di-<i>tert</i>-butylpyridine CAS RN: 585-48-8</p>
<p>D4652 200mg 1g</p>  <p><i>trans</i>-2,6-Diisopropyl-<i>N</i>-(2-pyridylmethylene)aniline CAS RN: 908294-68-8</p>	<p>B0468 25g 100g 500g</p>  <p>2,2'-Bipyridyl CAS RN: 366-18-7</p>	<p>D3134 1g 5g</p>  <p>4,4'-Di-<i>tert</i>-butyl-2,2'-bipyridyl CAS RN: 72914-19-3</p>	<p>P0221 1g 25g</p>  <p>1,10-Phenanthroline Monohydrate CAS RN: 5144-89-8</p>	<p>D0771 1g</p>  <p>Neocuproine Hemihydrate CAS RN: 34302-69-7</p>
<p>D0905 1g 5g</p>  <p>Bathophenanthroline CAS RN: 1662-01-7</p>	<p>T0361 25mL 100mL 500mL</p>  <p>Tributylphosphine CAS RN: 998-40-3</p>	<p>T2584 1g 5g</p>  <p>Tri-<i>tert</i>-butylphosphonium Tetrafluoroborate CAS RN: 131274-22-1</p>	<p>D4731 1g 5g</p>  <p>Di-<i>tert</i>-butyl(methyl)phosphonium Tetrafluoroborate CAS RN: 479094-62-7</p>	<p>T1165 25mL</p>  <p>Tricyclohexylphosphine (contains Tricyclohexylphosphine Oxide) (ca. 18% in Toluene, ca. 0.60mol/L) CAS RN: 2622-14-2</p>
<p>T2585 1g 5g</p>  <p>Tricyclohexylphosphonium Tetrafluoroborate CAS RN: 58656-04-5</p>	<p>D2411 1g 5g</p>  <p>Dicyclohexylphenylphosphine CAS RN: 6476-37-5</p>	<p>T0519 25g 100g 500g</p>  <p>Triphenylphosphine CAS RN: 603-35-0</p>	<p>T1024 5g 25g</p>  <p>Tri(<i>o</i>-tolyl)phosphine CAS RN: 6163-58-2</p>	<p>T2900 5g</p>  <p>Tris(4-fluorophenyl)phosphine CAS RN: 18437-78-0</p>
<p>T2526 1g 5g</p>  <p>Tris[3,5-bis(trifluoromethyl)phenyl]phosphine CAS RN: 175136-62-6</p>	<p>T1614 5g 25g</p>  <p>Tris(2,6-dimethoxyphenyl)phosphine CAS RN: 85417-41-0</p>	<p>T1643 1g 5g</p>  <p>Tri(2-furyl)phosphine CAS RN: 5518-52-5</p>	<p>T1666 1g 5g</p>  <p>Tri(2-thienyl)phosphine CAS RN: 24171-89-9</p>	<p>B1174 1g</p>  <p>1,2-Bis(dimethylphosphino)ethane CAS RN: 23936-60-9</p>
<p>B1137 10g 25g</p>  <p>1,2-Bis(diphenylphosphino)ethane CAS RN: 1663-45-2</p>	<p>B1138 5g 25g</p>  <p>1,3-Bis(diphenylphosphino)propane CAS RN: 6737-42-4</p>	<p>B1246 5g 25g</p>  <p>1,4-Bis(diphenylphosphino)butane CAS RN: 7688-25-7</p>	<p>B1960 1g</p>  <p>1,5-Bis(diphenylphosphino)pentane CAS RN: 27721-02-4</p>	<p>B1959 1g 5g</p>  <p>1,6-Bis(diphenylphosphino)hexane CAS RN: 19845-69-3</p>
<p>B3372 1g 5g</p>  <p>1,2-Bis(diphenylphosphino)benzene CAS RN: 13991-08-7</p>	<p>B2710 100mg 1g</p>  <p>1,1'-Bis(diisopropylphosphino)ferrocene CAS RN: 97239-80-0</p>	<p>B2711 100mg 1g</p>  <p>1,1'-Bis(di-<i>tert</i>-butylphosphino)ferrocene CAS RN: 84680-95-5</p>	<p>B2027 1g 5g 25g</p>  <p>1,1'-Bis(diphenylphosphino)ferrocene CAS RN: 12150-46-8</p>	<p>B2383 5g 25g</p>  <p>(+/-)-BINAP CAS RN: 98327-87-8</p>

B2867 5g 25g  DPEphos CAS RN: 166330-10-5	B2709 1g 5g 25g  Xantphos CAS RN: 161265-03-8	B5239 200mg 1g  4,5-Bis(dicyclohexylphosphino)-9,9-dimethylxanthene CAS RN: 940934-47-4	B2717 100mg 1g  4,6-Bis(diphenylphosphino)phenoxazine CAS RN: 261733-18-0	D3389 1g 5g  DavePhos CAS RN: 213697-53-1	
B2630 100mg 1g  2,2'-Bis(diphenylphosphino)biphenyl CAS RN: 84783-64-2	B5957 100mg 500mg  2,2'-Bis[bis(3,5-dimethylphenyl)phosphino]-1,1'-biphenyl CAS RN: 325773-62-4	B4595 1g  1,3-Bis[(di-tert-butylphosphino)oxy]benzene CAS RN: 338800-20-7			
<h2>Directing Group Introducing Agents</h2>			A0419 5g 25g 100g  8-Aminoquinoline CAS RN: 578-66-5	D4264 1g 5g  2-(Diisopropylsilyl)pyridine CAS RN: 1232692-92-0	P1902 1g  2-(1H-Pyrazol-5-yl)aniline CAS RN: 111562-32-4
<h2>Additives</h2>			B0089 25g 100g 500g  p-Benzoquinone CAS RN: 106-51-4	T1244 25g 500g  p-Toluquinone CAS RN: 553-97-9	D0686 1g 5g 25g  p-Xylquinone CAS RN: 137-18-8
D2234 1g 5g 25g  2,6-Dimethyl-1,4-benzoquinone CAS RN: 527-61-7	D3411 100mL  Di-tert-butyl Peroxide CAS RN: 110-05-4	B3153 100g  tert-Butyl Hydroperoxide (70% in Water) CAS RN: 75-91-2	T1560 5g 25g  TEMPO Free Radical CAS RN: 2564-83-2	F0335 5g 25g  NFSI CAS RN: 133745-75-2	
P1015 5g 25g  Koser Reagent CAS RN: 27126-76-7	I0330 10g 25g 250g  Iodobenzene Diacetate CAS RN: 3240-34-4	D0798 25g 500g  Dimethyl Sulfoxide CAS RN: 67-68-5	T0431 25g 100g 500g  Trifluoroacetic Acid CAS RN: 76-05-1	O0310 25g 500g 2KHSO₅ · KHSO₄ · K₂SO₄ Potassium Peroxymonosulfate CAS RN: 37222-66-5	
L0224 25g 500g  Lithium Carbonate CAS RN: 554-13-2	S0560 300g  Sodium Carbonate CAS RN: 497-19-8	P1748 300g  Potassium Carbonate CAS RN: 584-08-7	C2160 25g 100g  Cesium Carbonate CAS RN: 534-17-8	T2052 100mL 500mL TiCl₄ Titanium(IV) Chloride (14% in Dichloromethane, ca. 1.0mol/L) CAS RN: 7550-45-0	
T3238 100mL 500mL TiCl₄ Titanium(IV) Chloride (ca. 19% in Toluene, ca. 1.0mol/L) CAS RN: 7550-45-0	S0463 5g 25g AgSbF₆ Silver Hexafluoroantimonate(V) CAS RN: 26042-64-8	S0898 1g 5g Ag⁺ (CF₃SO₂)₂N⁻ Silver Triflimide CAS RN: 189114-61-2	S0978 5g 25g  Sodium Pivalate Hydrate CAS RN: 143174-36-1	P2354 5g 25g  Potassium Pivalate CAS RN: 19455-23-3	



Ordering and Customer Service

TCI AMERICA

Tel : 800-423-8616 / 503-283-1681
Fax : 888-520-1075 / 503-283-1987
E-mail : Sales-US@TCIchemicals.com

TCI EUROPE N.V.

Tel : +32 (0)3 735 07 00
Fax : +32 (0)3 735 07 01
E-mail : Sales-EU@TCIchemicals.com

TCI Deutschland GmbH

Tel : +49 (0)6196 64053-00
Fax : +49 (0)6196 64053-01
E-mail : Sales-DE@TCIchemicals.com

Tokyo Chemical Industry UK Ltd.

Tel : +44 (0)1865 784560
E-mail : Sales-UK@TCIchemicals.com

TCI Chemicals (India) Pvt. Ltd.

Tel : 1800 425 7889 / 044-2262 0909
Fax : 044-2262 8902
E-mail : Sales-IN@TCIchemicals.com

梯希爱(上海)化成工业发展有限公司

Tel : 800-988-0390 / 021-67121386
Fax : 021-6712-1385
E-mail : Sales-CN@TCIchemicals.com

TOKYO CHEMICAL INDUSTRY CO., LTD.

Tel : +81 (0)3-5640-8878
E-mail : globalbusiness@TCIchemicals.com

Availability, price or specification of the listed products are subject to change without prior notice. Reproduction forbidden without the prior written consent of Tokyo Chemical Industry Co., Ltd.